

# How to install and activate the MestreNova NMR Campus licence

## Introduction

The NMR campus licence, of MestreNova software, can be installed and activated from any devices providing that you are using the university network (wired network, Wi-Fi or VPN). Licence is perpetual but requires to connect to the university network every 90 days to remain in active status.

The actual campus NMR licence is inclusive of:

- NMR
- qNMR (quantitative NMR)
- RM (reaction monitoring for array and kinetics experiments)

## Licences file


The campus NMR licence is sent by e-mail in zip format.

## How to instal the program


1. Go to the MestreLab internet page <https://mestrelab.com/download/mnova/> and download the MestreNova version suitable for your operative system. DO NOT download the last version available, but the 15.0.0 version;
2. Launch the installation file and a window will appear: go on without customize the installation. At the end of the installation, click on "Finish" button as in the picture below;




3. Launch the program: a window asking for the licences will appear (see picture below):

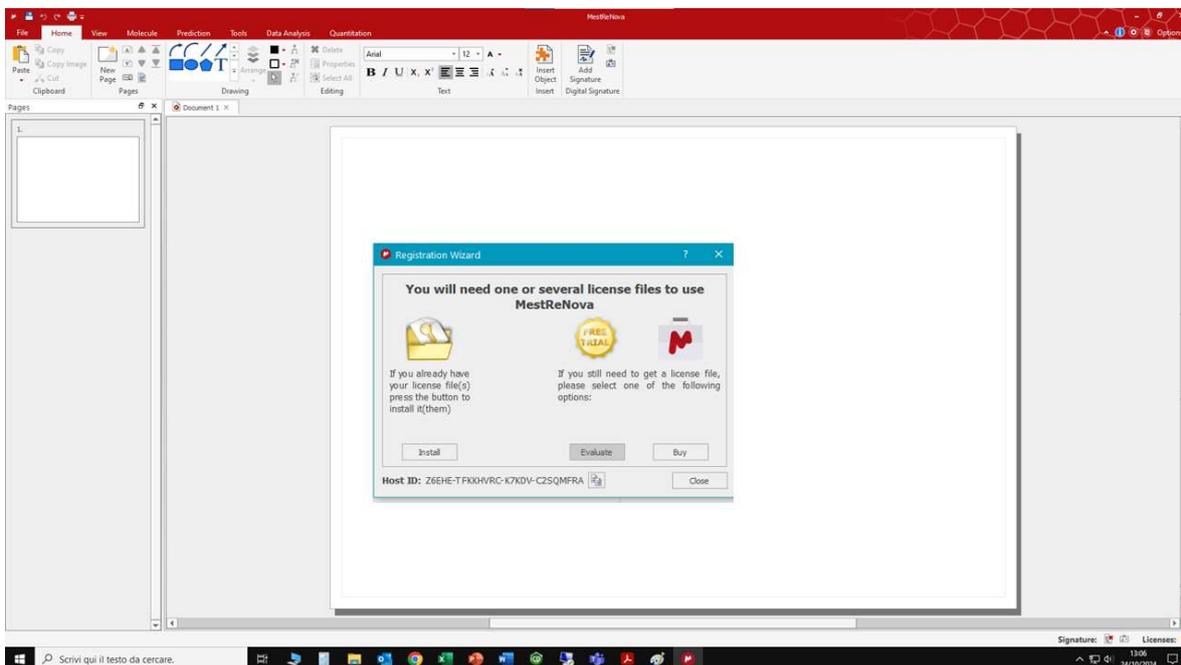

4. Click on the “Install” button and select the path where you have saved the licences and install them one by one (in the folder there are 3 licences: it possible to install all of them or only the ones of interest);

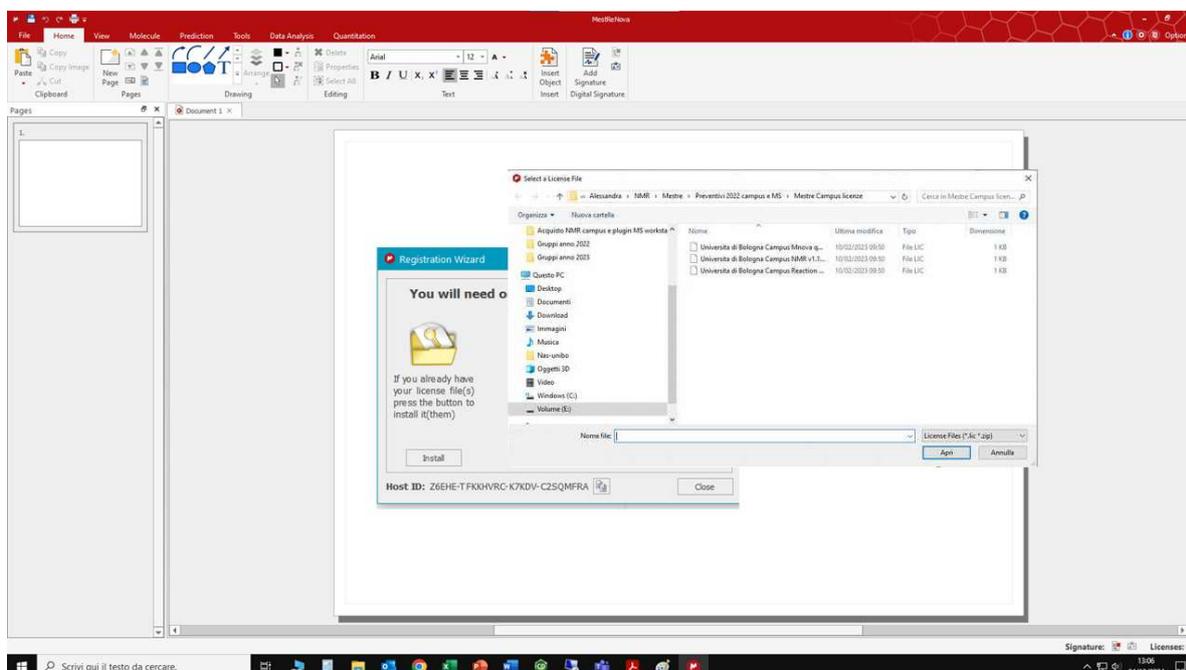

5. Close and restart the program. To verify if the licences are correctly installed, click on the second to last icon on the red bar on the right-hand side and see if all licences are activated as in the picture below;

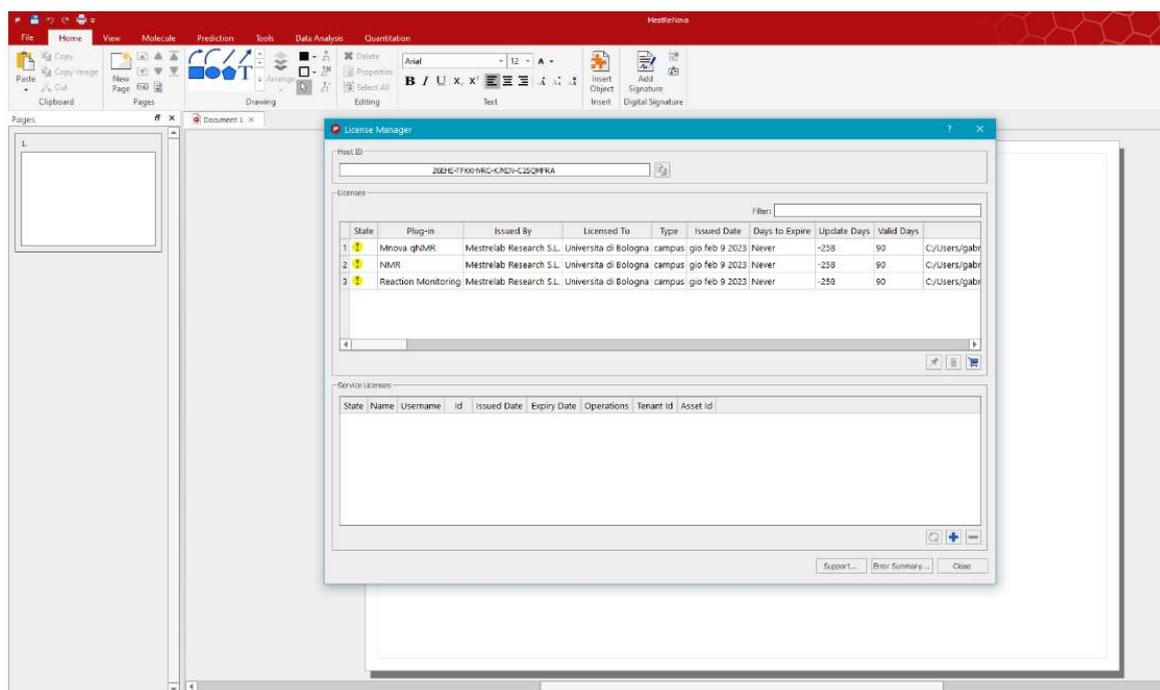

**P.S.:** The presence of the 3 exclamation points does not indicate a problem with the licences, but only that the licenses are not up to date.